



Bios

Driving a Higher Standard
in Flow MeasurementSM

Bios Optimizer Collect Upgrade Procedure

To upgrade to Bios Optimizer Collect from Bios Optimizer Collect light requires:

- The purchase of a CD key from Bios for Optimizer Collect, price \$495
- Bios will e-mailed you the CD key
- Start Collect light on your PC
- On the Collect light start screen, click on Registration>Bios Optimizer Collect light
- Record the Registration Code
- Go to the Bios Optimizer Collect activation page and enter the required information- <http://www.biosint.com/Products/OptRegistration.php>
- You will be sent a Activation Code via e-mail
- On the Collect light page, click on Registration>Bios Optimizer Collect light and enter your new Activation Code